


Applying a Marker Bar

1. Select **Measure/ Calibration/ Select Spatial** -This opens the below dialog – select the relevant calibration

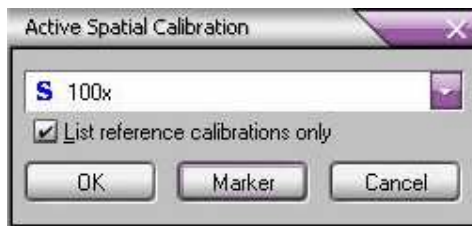

2. Select **Marker** – this opens the Marker Dialog box



3. **Enter a Marker Bar length** – our suggestions are below

Objective Power	Suggest Marker Size
4x	250 microns
10x	100 microns
20x	50 microns
40x	25 microns
60x	20 microns
100x	10 microns

4. **Select OK** – the marker will now be placed on your image using the left mouse button move it to the correct position and the right click when happy.

The marker bar can now be saved along with your image