

Opening Images as a Sequence

1. Select **File/Open** to open the regular **Open File** dialog
2. At the bottom right of the dialog select **Open as Sequence**
3. Select the last image in the sequence then select **“Shift”** and highlight all other images you wish to open
4. Select **Open**

Your image will now be open as a sequence file and can be saved as either a .seq or .avi file.

Alternative Method

This is more common in older versions of Image-Pro (pre version 6.0)

1. Select **Sequence/ Merge Files** – this will open the Open File Dialog
2. Select the last image in the sequence then select **“Shift”** and highlight all other images you wish to open
3. Select **Open**

